

	<p>blue-red colonies with gas regardless of size or intensity of colour or red colonies with gas. Typical E.coli colony morphology is blue to blue-red colonies with gas regardless of size or intensity of colour. Decimal dilutions of the poultry or poultry products can also be prepared and tested.</p> <p>A count of colonies on each plate is then performed, the result is then calculated taken into account the dilution factor and poultry portion/mass. The number of micro-organisms per g or cm² is then calculated.</p>
Reporting Unit	E.coli Count cfu/g, cfu/bird or cfu/cm ²
LOR/LOQ	< 10

Minimum amount of sample required for analysis	Poultry unit	Turnaround time	3 days
--	--------------	-----------------	--------

Author:	Document Controller	Date:	19 th December 2017
Authorised By:	National Quality Manager	Date	19 th December 2017

CONFIDENTIAL